

NMR data access at **imsercdata.northwestern.edu** (*archived and real-time data*):

To view your group's archived data, you must complete the training at:

<http://129.105.219.102/GettingStarted/multiscreen.html> and agree to follow IMSERC's policies. It typically takes less than one week to activate an account. If you can't login with your account after it is activated, try changing your NUID password at: <https://validate.it.northwestern.edu/idm/user/login.jsp>

For windows users:

At "Map network drive" dialog box, input `\\imsercdata.northwestern.edu\share`, use your netid and passwd or ad\netid and passwd when promoted. "share" is your PI's last name, all lower case (see last paragraph for complete list).

For mac users:

"Apple+k" to open "connect to server" dialog box, input `smb://imsercdata.northwestern.edu/share`, the rest is same as windows.

All IMSERC data is organized into five subdirectories (ms, nmr, nmr_rt, xrd, others). Click "nmr" for archived and "nmr_rt" for real-time nmr data from all NMR instruments for your group.

NMR data access with sftp at **wakko3** (*real-time data*)

Get winscp at <http://winscp.net/eng/download.php>
or filezilla at <https://filezilla-project.org/download.php>

Setup the program to access **wakko3.chem.northwestern.edu** (real time access) with *sftp* protocol. Once you are connected to wakko3 (user: walkon, passwd: GO*****), all the nmr instruments are under /nfs. You can also get full set of vnmrj manuals at /vnmr/help/pdf for your reference if you want.

Available shares to map:

Everyone has access to: public, scratch

You must request access to: ameer beddows cmidd dravid freedman gaillard geiger gray grzybowski harris hersam high_throughput ho hodgson hoffman hupp hurtgen ibers jewett joester kanatzidis kelleher kung lewis marks meade messersmith mirkin mondragon mrksich nana nguyen northrup notestein ohalloran ohmx poeppelmeier priest radhakrishnan rosenzweig scheidt shull silverman stair statsyuk stoddart stupp thomson torkelson tyo underedu vanduyne wasielewski watterson weiss weitz